

coolingtech

Digital Microscope User Manual

English

Thank you for purchasing our Digital Microscope

It is a slim type tool and can be easily used in different fields.
Including:

1. Textile industrial for textile inspection
2. Printing inspection
3. Industrial inspection: PCB, Precision machinery...
4. Educational purpose
5. Hair examination
6. Skin examination
7. Microbiological observation
8. Jewelry & coin (Collections) inspection
9. Visual Assistance
10. Others

Currently we have two models of Digital Microscope: USB port and TV port formats.

The USB microscope must be connected to the computer or mobile phone, tablet PC. and TV microscope be connected to visual device (N/P). User interface is very friendly and you can do the Measure work easily.

You can find more information of the product on this manual.

CONTENTS

Introduction	2
Parts of Microscope	3
Specification	4
Product Outlook & Standard parts.....	4
Notes before use.....	4
Hardware and system requirements	4
How to install the Driver software	5
How to install the measurement software	10
Function options	10
How to do the calibration	14

New features of Coolingtech software version 4.5 :

- All functions support win10 system
- Software supports 19 languages (Chinese, English, Korean, Japanese, Spanish, German, French, Russian, Thai, Italian, polish, Portuguese, Dutch, Turkish, Malay, Greek, Persian, Arabic)
- Screenshot interface sort by chronological order from the back to the front

Parts of Microscope



Specification

Sensor: high performance chip
Main control chip: dedicated master 24Bit DSP
Photo / video: software control
Light source: 8 LED white lights
Static resolution: SD 640*480; HD 1200*1600; ultra clear 2592*1944
Digital zoom: 5X
Working distance: manual adjustment 0~40mm
Image resolution: SD 640*480P, HD 1280 x 720P; Full HD 1960*1080P
CD:driver software, measurement software, use manual
Support operating system: 32bit and 64bit Windows,, XP, VISTA, WIN7, WIN 8, WIN 10; MAC, Linux

Power supply: USB (5V DC), WIFI or mobile OTG cable

Interface: USB 3.0 & USB 2.0 & USB 1.1 compatible, OTG

Dynamic Frames: 30f/s Under 600 LUX Brightness

Illumination range: 0 ~ 30000LUX adjustable by USB cable Hardware requirements: Unlimited

Language support: [Chinese \(Simplified\)](#), [Chinese \(traditional\)](#), [English](#), [Korean](#), [Japanese](#), [Spanish](#), [German](#), [French](#), [Russian](#), [Thai](#), [Italian](#), [polish](#), [Portuguese](#), [Dutch](#), [Turkish](#), [Malay](#), Greek, Persian, Arabic (other languages need to customize)

Product color: matte black, other colors can be customized

Body size: 112 mm (length) 33 mm (outer diameter)

Single package weight: 265g

product support [Windows 2000](#), [XP](#), [VISTA](#), [WIN 7](#), [WIN 8](#), [WIN 10](#) Pentium 1G, Celeron, AMD 1G & above, 128MB Memory, 150MB Hard Disc memory space, 16-bit & above VGA, CD-ROM, USB2.0 or USB1.1

Product Outlook & Standard parts

1. Digital Microscope (1pcs)
2. USB 2.0 cable
3. Metal fixed Stand (1pcs)
4. CD (Driver software, Measurement software, User Manual)
5. Microscope Micrometer Calibration Rule


Notes before use

1. Don't disassemble the digital Microscope or change the interior parts, it can cause damage.
2. Don't clean the Microscope with alcohol organic solvents
3. Don't touch the lens with your fingers.
4. Avoid outdoor use if possible.
5. Storage temperature , 0°C ~ 40°C, Humidity: 45%RH ~ 85RH%.
6. In case the product gets wet, leave PC connection immediately. And do not disassemble or dry by hair dryer. Send to repair center if the digit microscope was effected by liquid or other elements.
7. Measurement Data only for reference

Hardware and system requirements

Digital Microscope is a free Driver Product, it can be worked immediately after connected to the computer with operation system above Win2000 version.

1. Connect the Portable digital microscope by USB cable to your computer, double click 'my computer' icon on your computer desktop, below interface will be shown up.

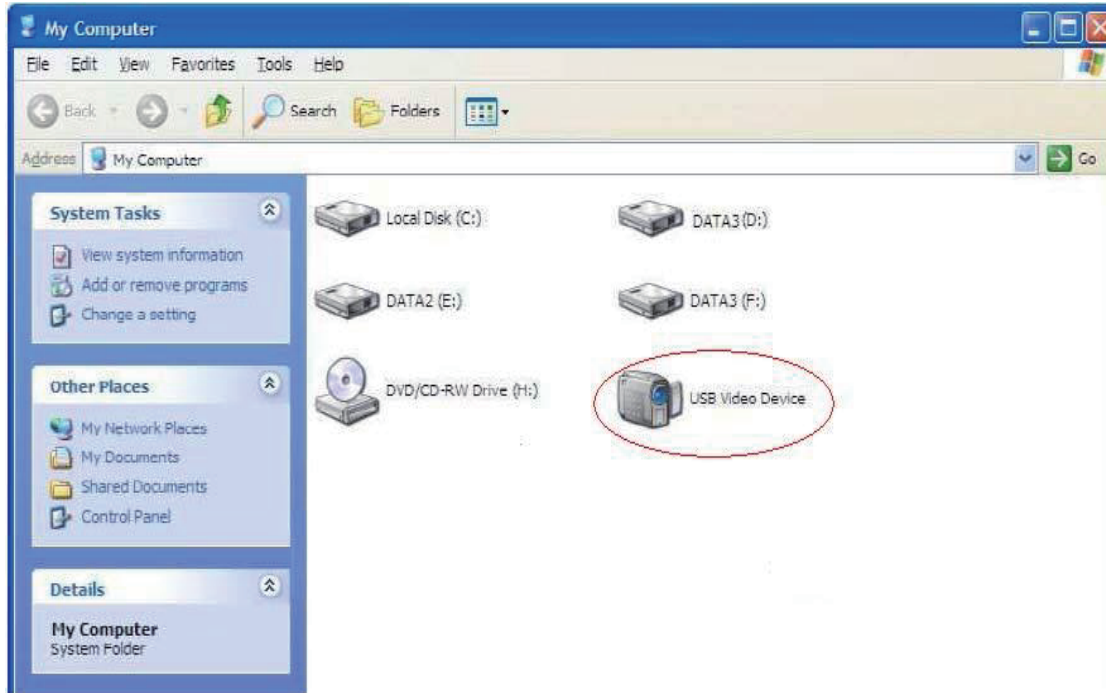

2. Double click the Video device icon as above red marked in your computer. Now it is ready to use.


If you can't find the Video device, or if your computer can't recognize it, please follow below steps to install the driver.

Insert the attached CD, system will auto run to the interface as below and follow the steps to complete the installation:
(NOTE: If the computer can't auto run the disc, you need to operate manually by click "My Computer" → "DVD/CD Driver" → run "AUTORUN.EXE")



How to install the Driver software


Click **Driver** to continue, the screen will show the installation progress.



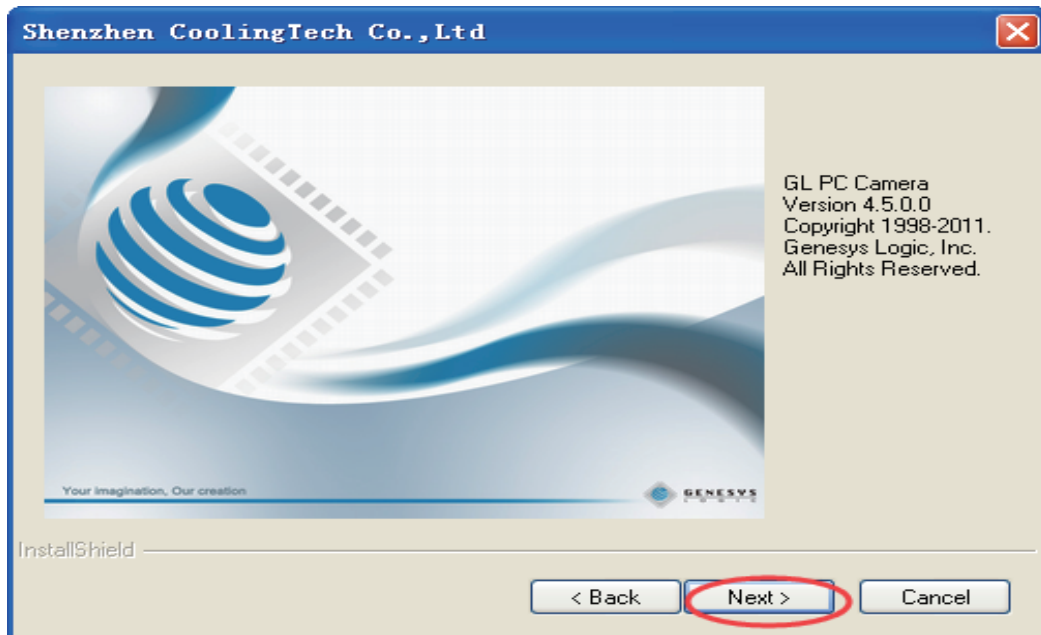


Click“Next”——Click“Next”——Click“Install”——Click“是(Y)”——Click“Finish”

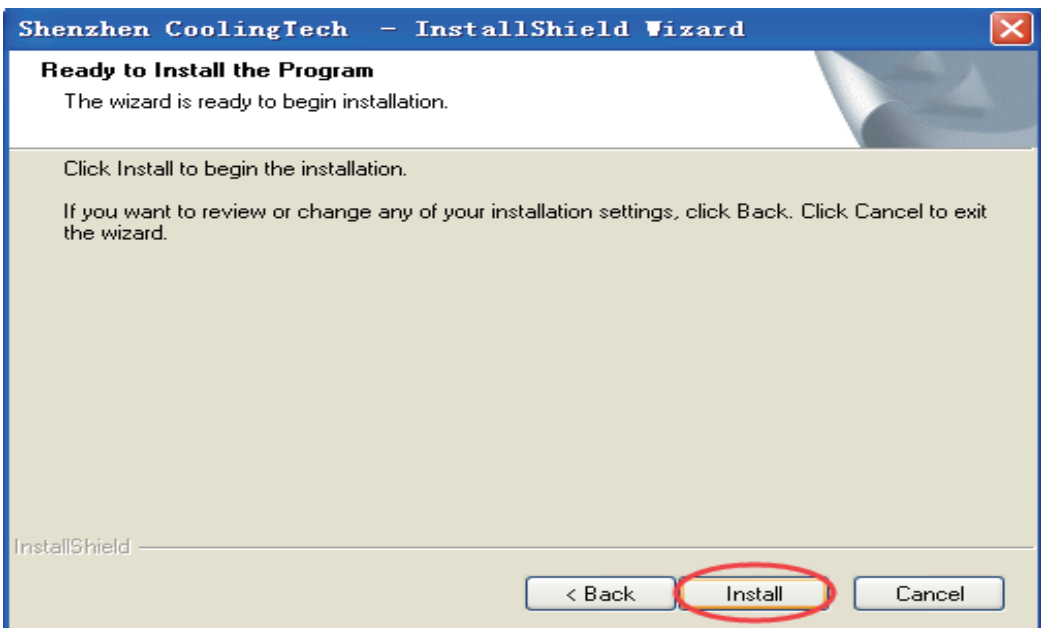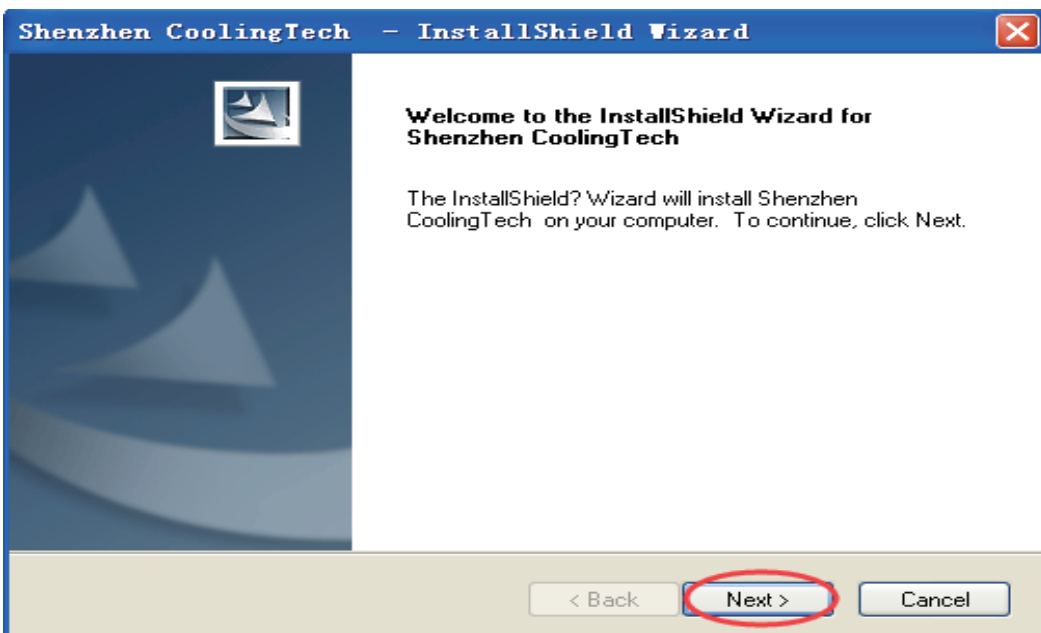

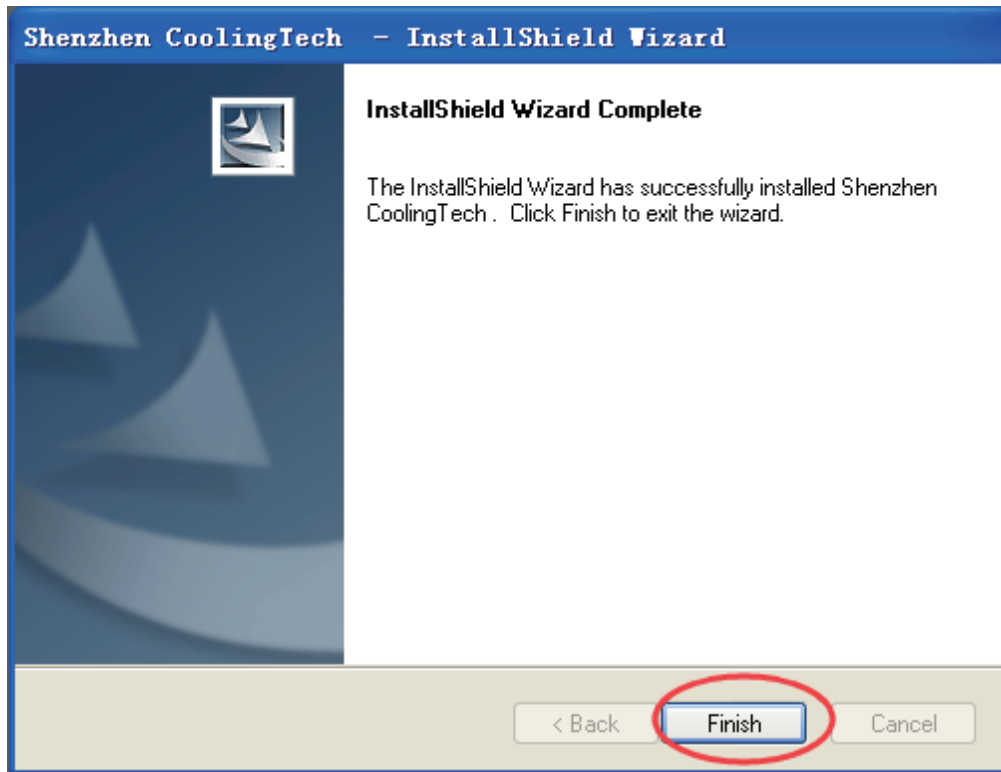


Select 'Finish' at last. Driver installation completed, if it is the first installation, it is recommended to restart your computer. If warning information appear, please click "continue".

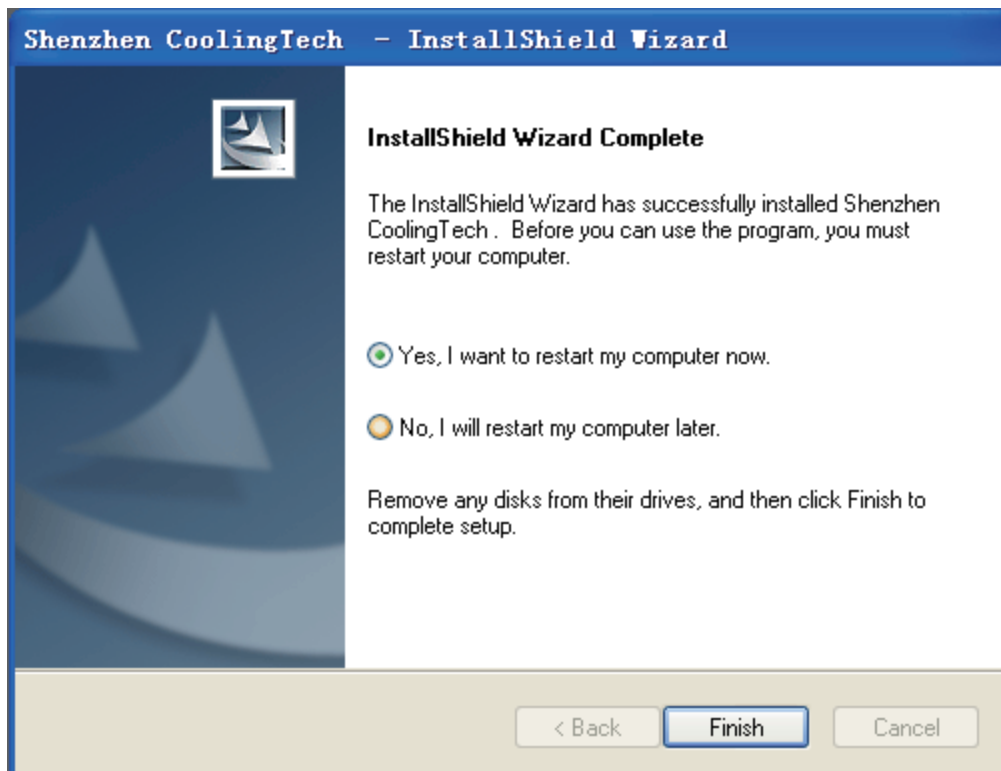




After the installation is complete, the following icon appears on the desktop, double-click the icon, click on the "device" option, select "USB 2 PC camera", will see the picture of the microscope capture

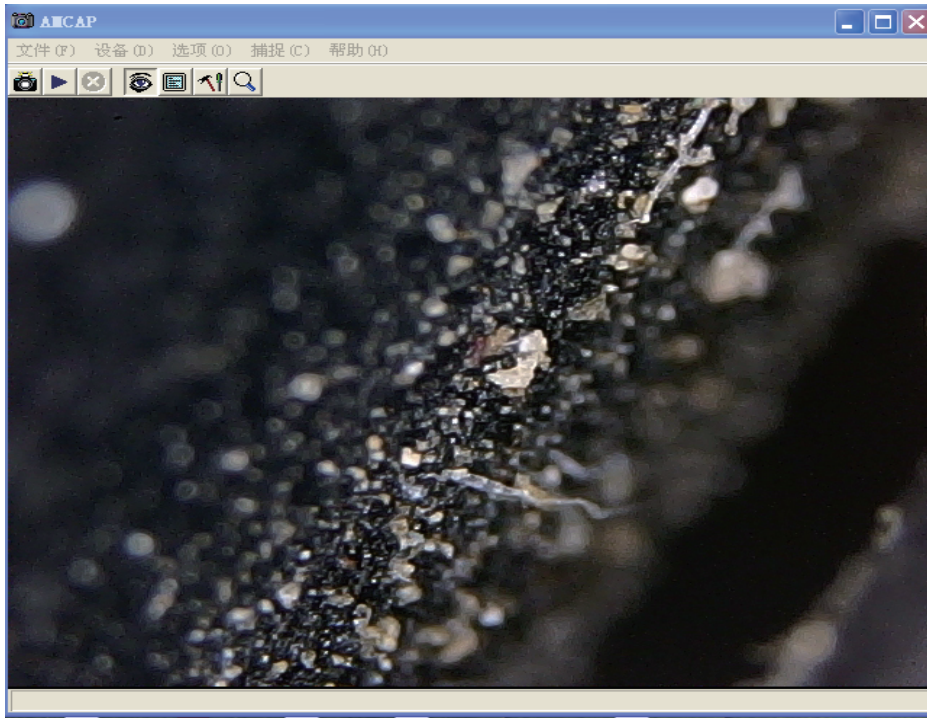

options in Menu bar can be set according to needs, for example, click the "options" - "set the video effect", in dialog box, you can adjust the picture effect. If you need to restore default settings, click "default".

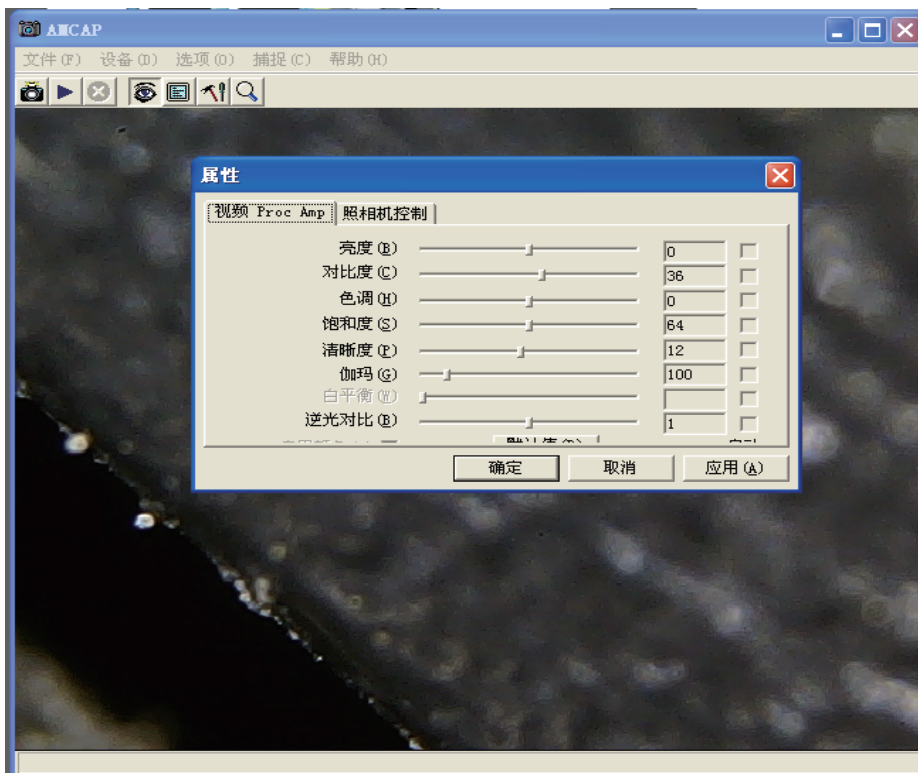




note: in the amcap software, the 'ZOOM' button can increase the magnification 5 times;and double click or long press 'SNAP' button can take photo.Set photo save path:secler 'capture', 'set image path'.
The amcap software can do general observation, if you need to measure objects,please install "Coolingtech" software in CD. installation steps as follows:

How to install the measurement software




Please connect the product to your computer before installation. click measurement **Measurement** , or open CD, double click Measurement.exe, select 'next' → setting setup path, select 'next' → select 'close'. In the installation process, if system prompt need to install 'Net framework V2.0', then open the 'dotnetfx.exe' in CD, to install Net framework V2.0, then you can successfully install measurement software.

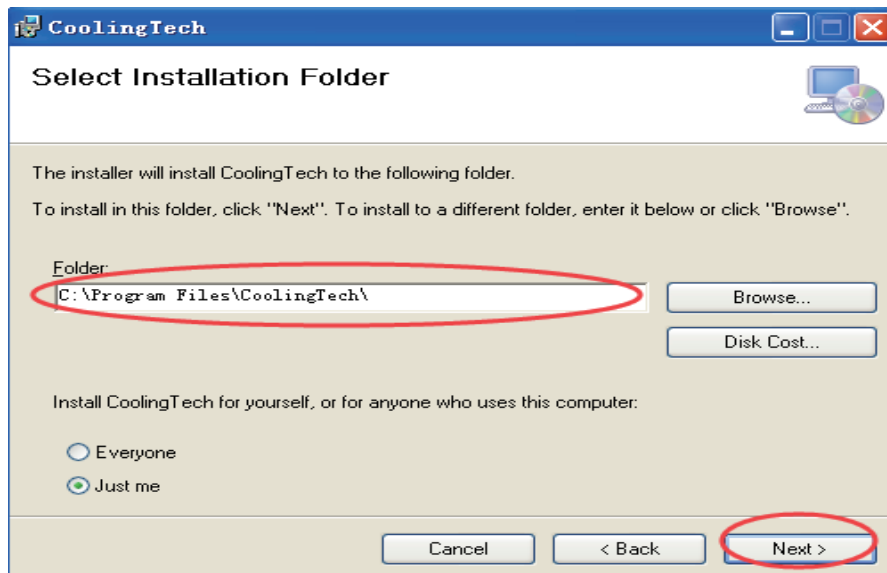

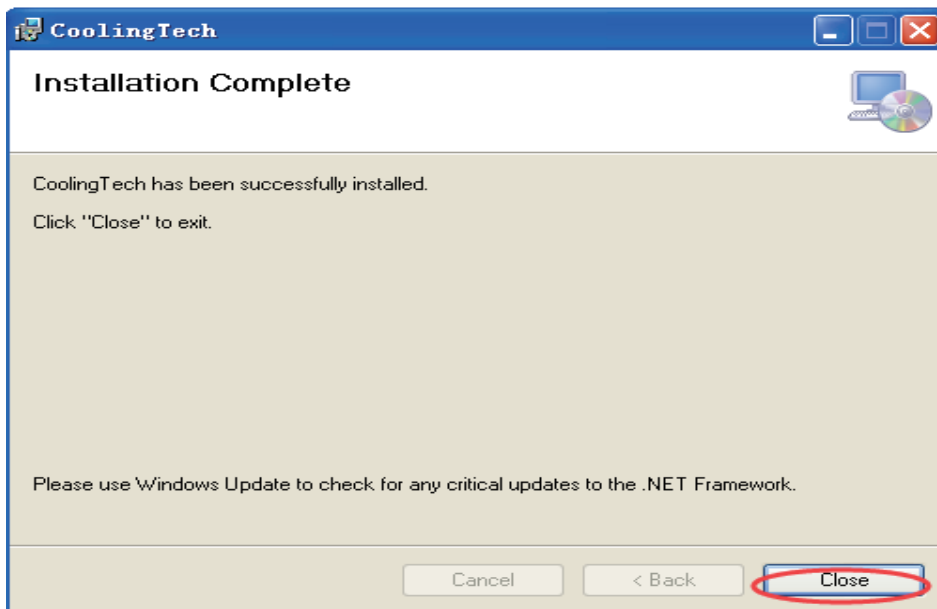
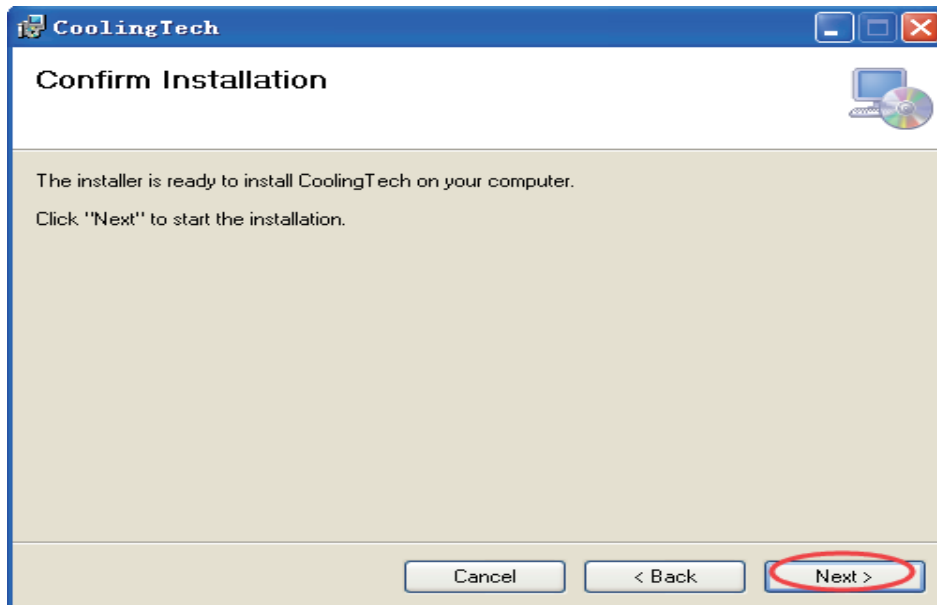


After the successful installation of the program, click on the desktop called "CoolingTech MicroScope" folder, enter the following interface. Double click the 'CoolingTech MicroScope' Icon.

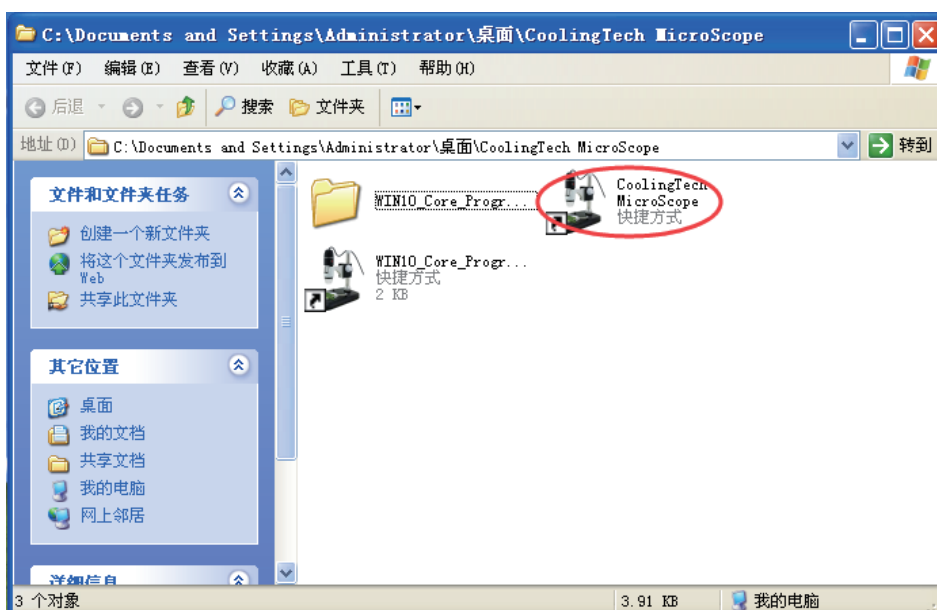





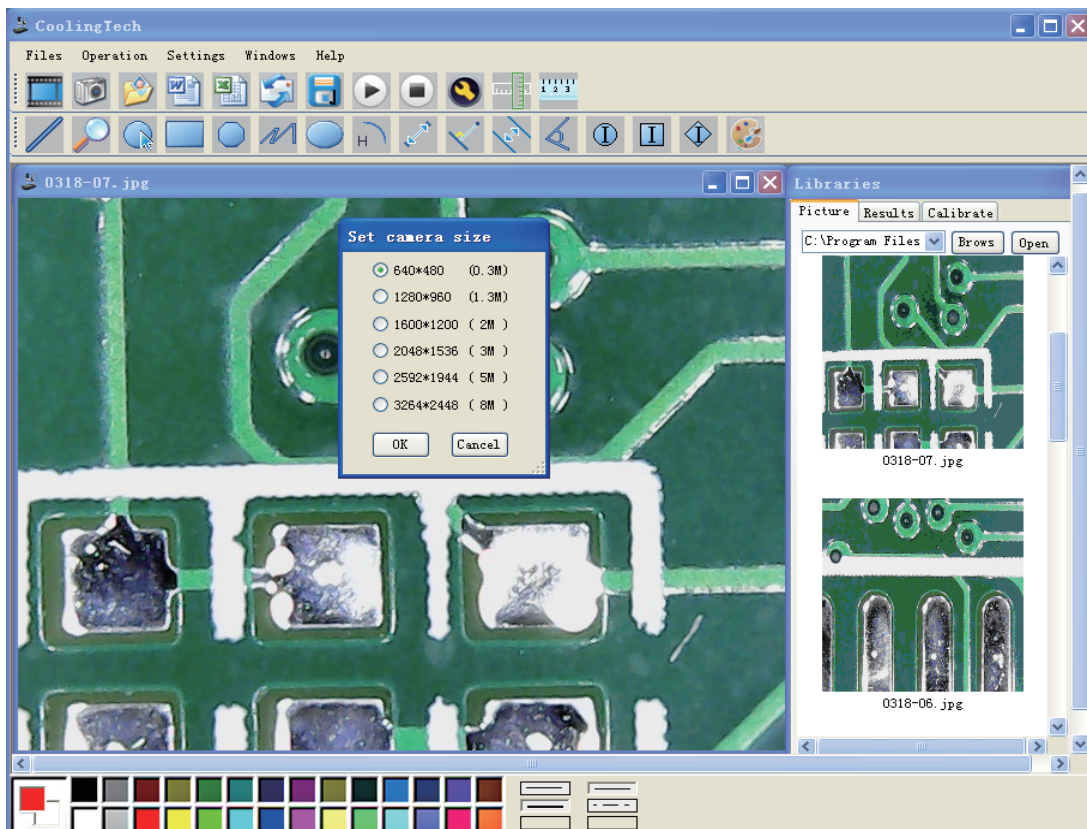
Follow the prompts to connect the device, the program automatically runs to the main interface.

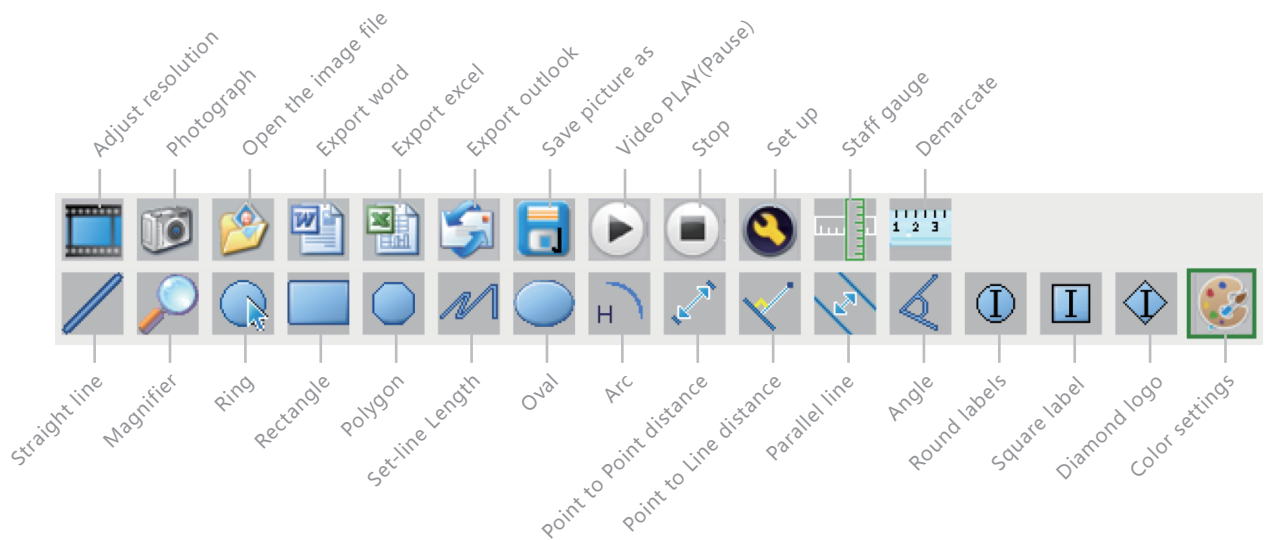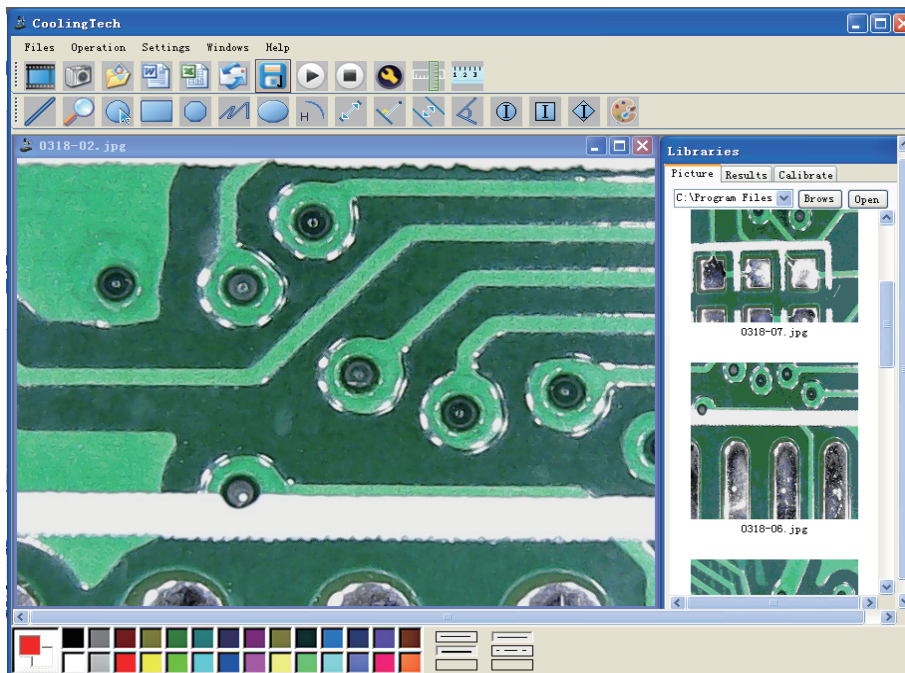


19 languages are optional: "set" - "language" - (select the language you want)

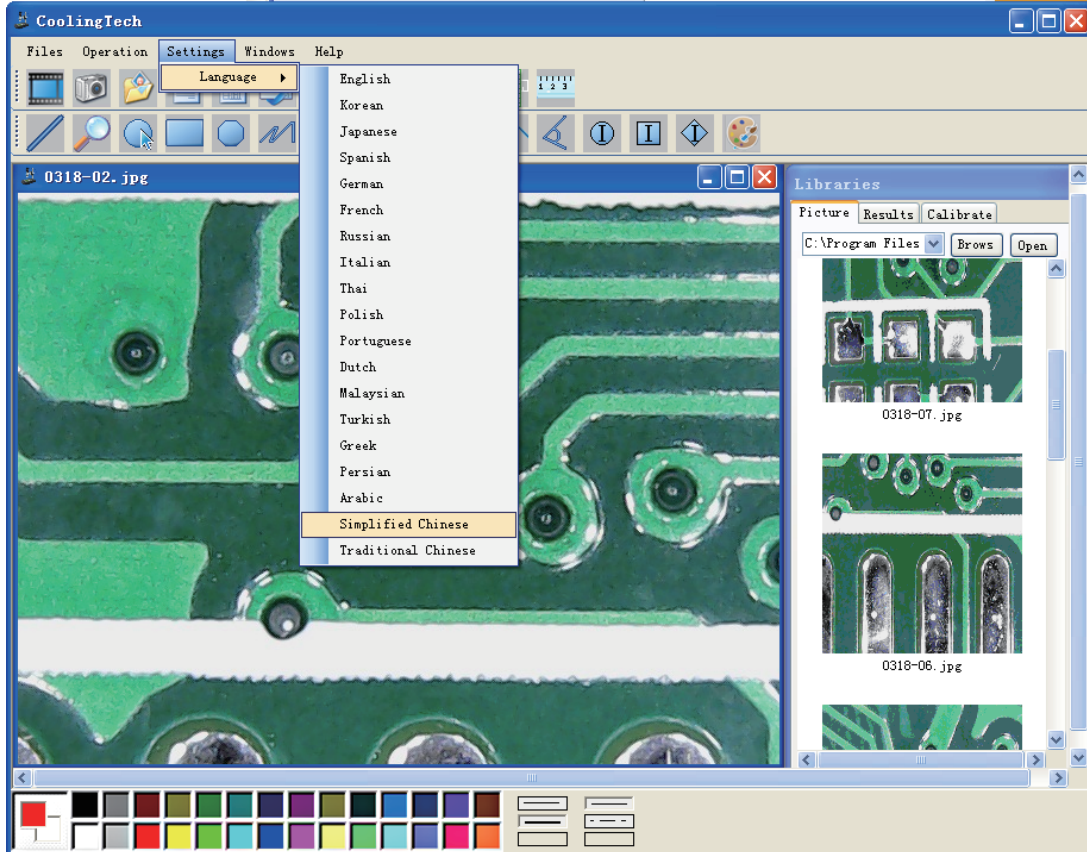


Click  , select the video resolution, video resolution range is 320*240 to 1960*1080, the default resolution is 640*480. click  to take pictures. The latest pictures at the top.




How to do the calibration

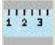
Put the microscope on the calibration ruler (gift with product) , adjust the roller to get a clear picture, press pause, click the icon  to do calibration. In order to improve the measurement accuracy, you must do calibration before using the microscope to measure. the magnification change each time, the calibration should be changed. It is recommended to measure the same actual value several times, and then take the average value.


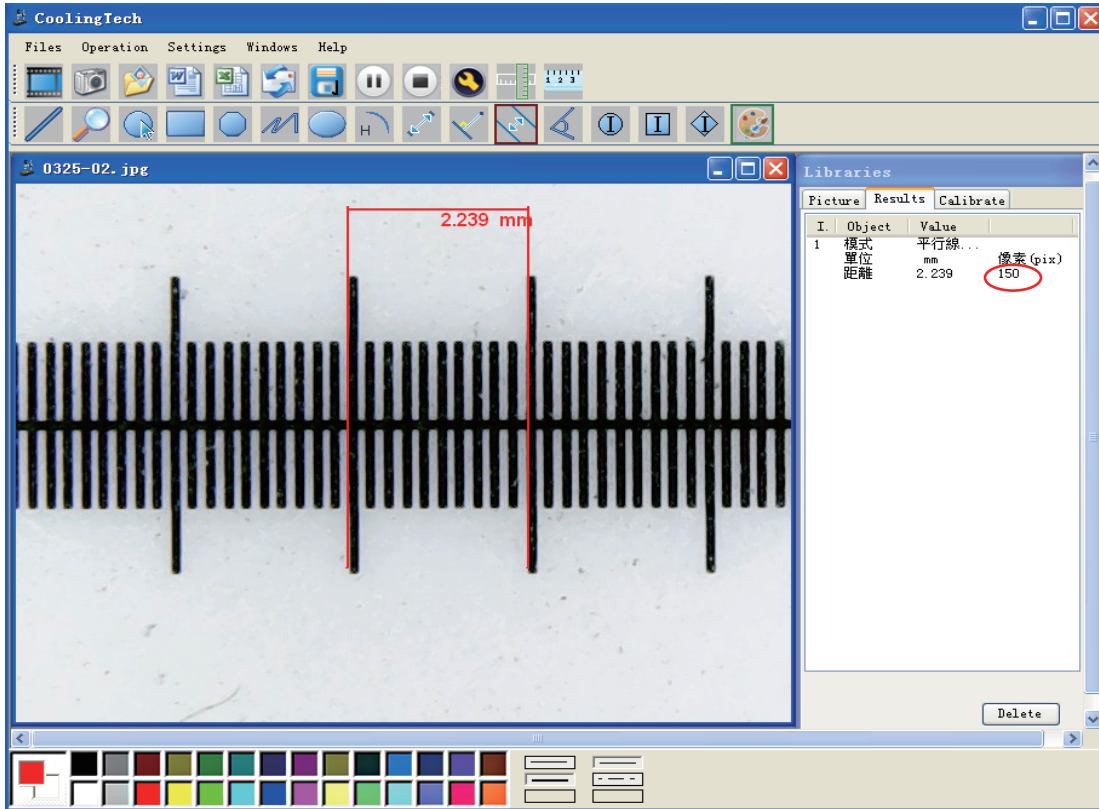


Enter the actual distance value, and the corresponding unit. The calibration is done. This calibration can be used to measure at this magnification.


Note: magnification change, must be corrected, calibration result will be different at the same time.

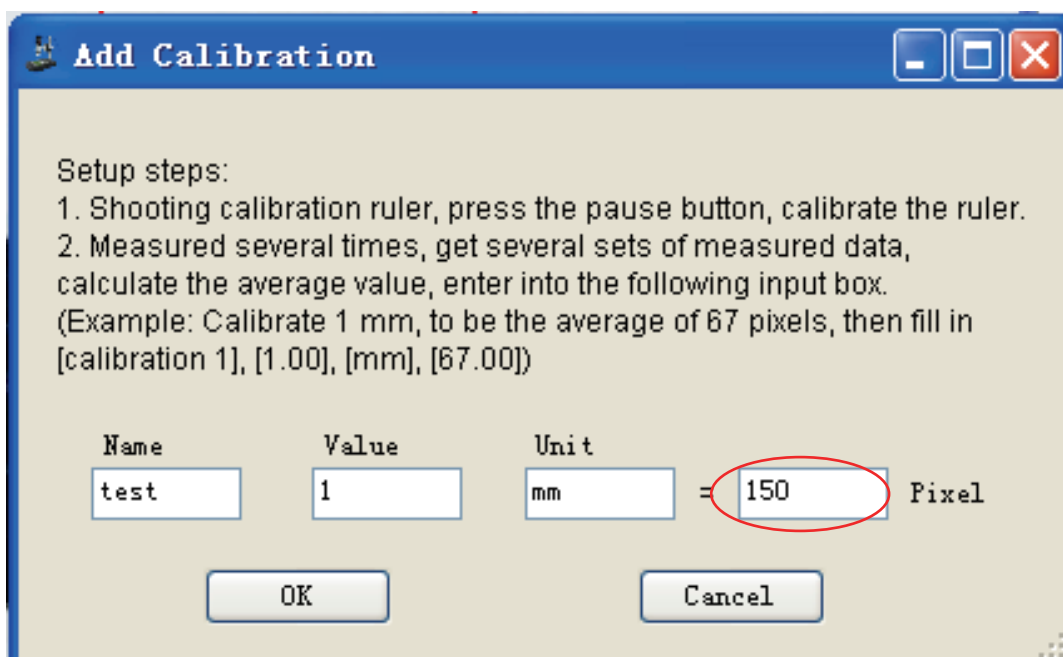


Calibration done, do not make any changes to the microscope, replace object with the calibration ruler and put it in the right position before the microscope. Click  to take pictures. Double click on the image, then you can measure it. The software can measure various points, lines, shapes, length, angle, radian and other parameters. Measurement results can be exported by Word or Excel, mail.

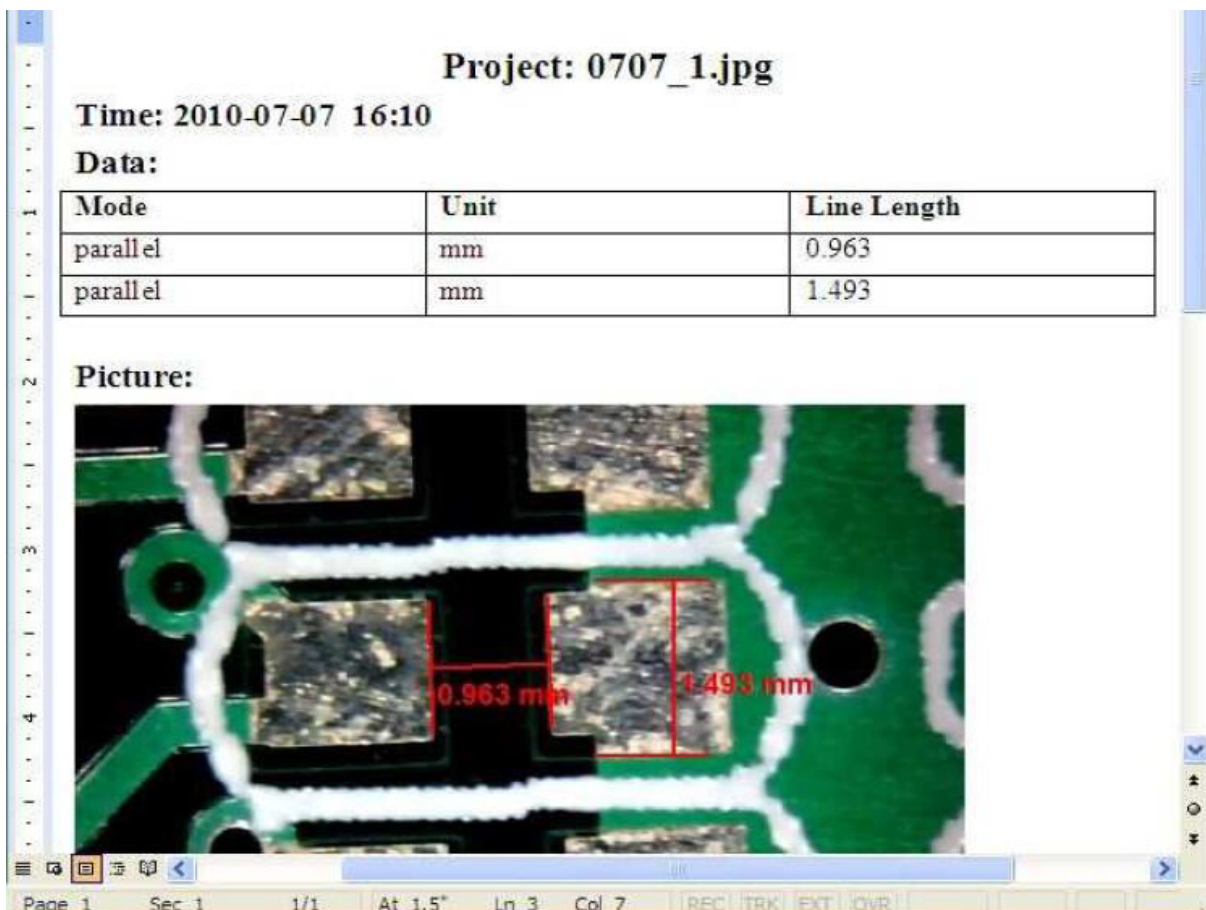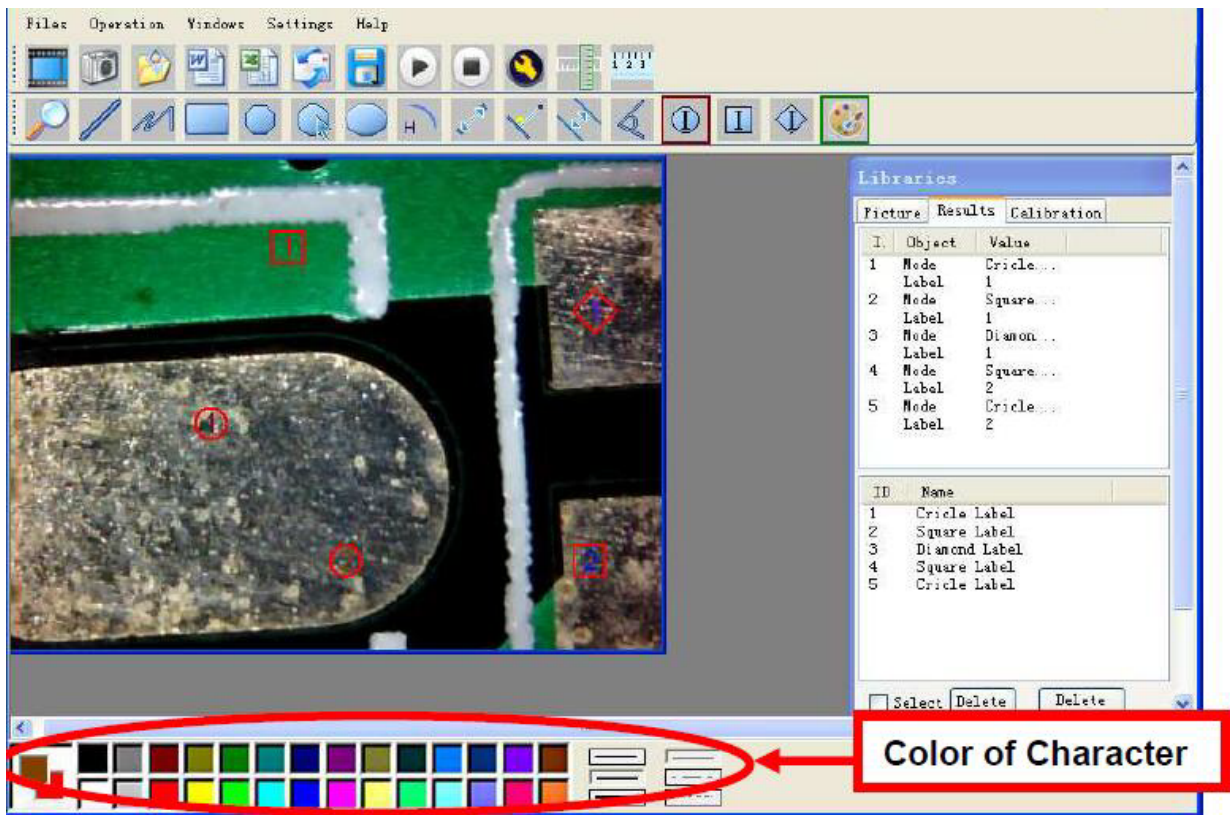

Click  , set screen brightness, contrast and other parameters.

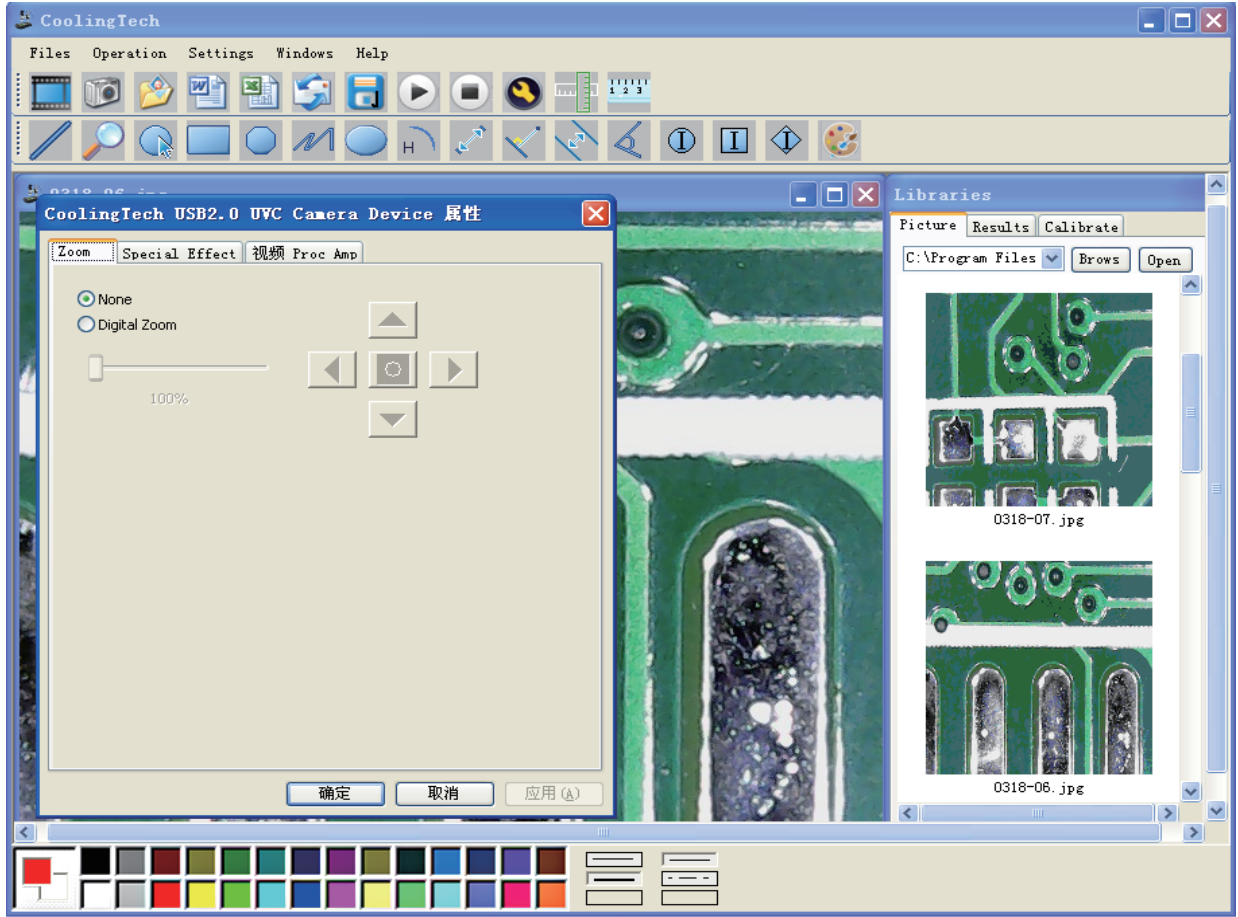


Click  Ruller, as followings

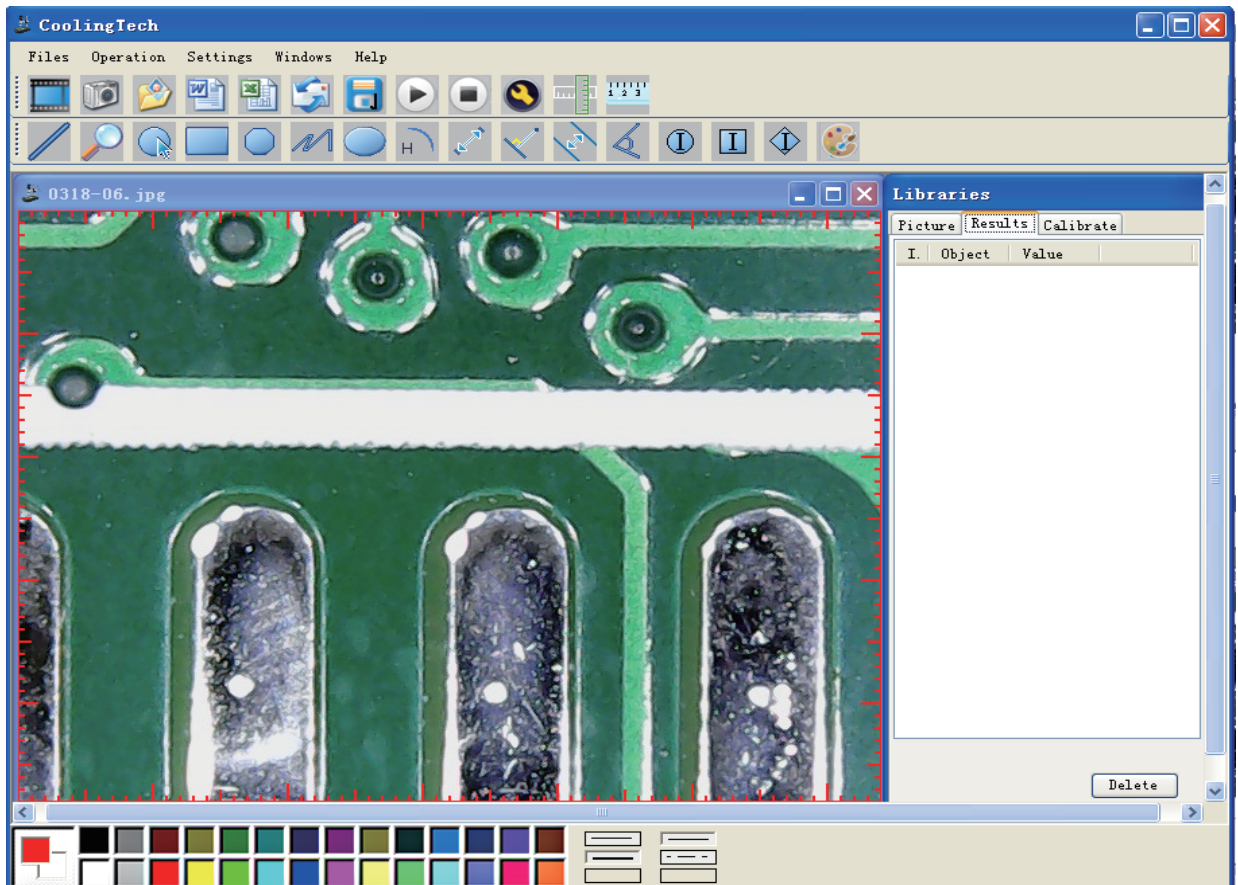




Annotation and color selection

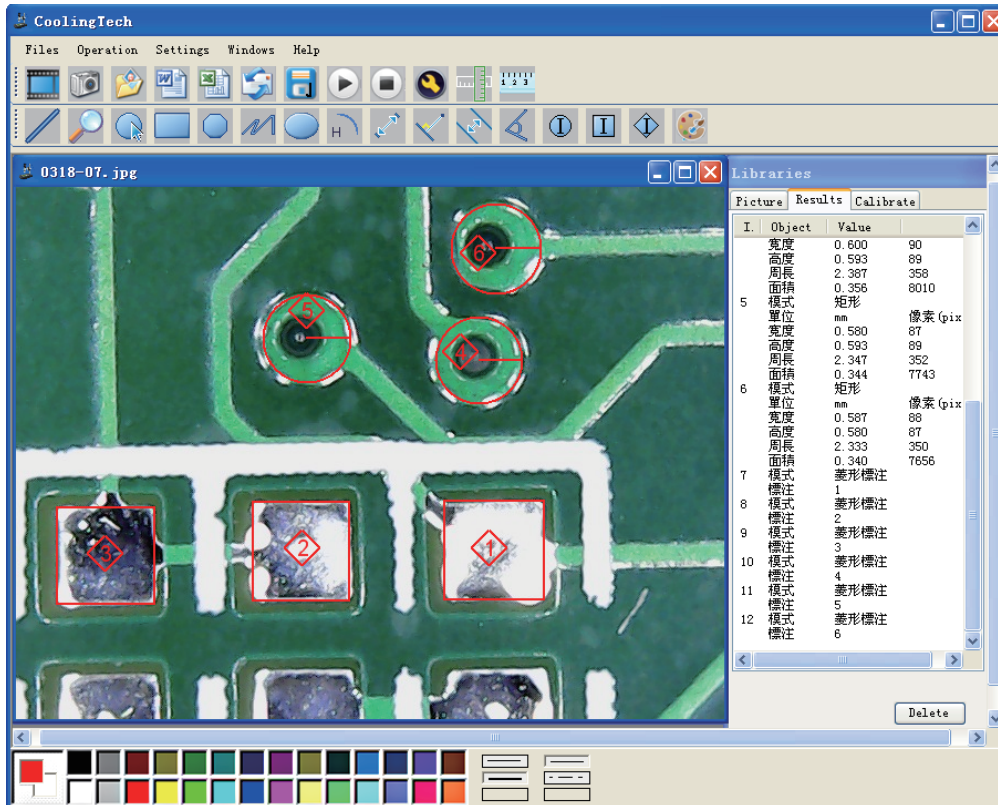


Icon description(left to right):


Linear measurement, circular measurement, rectangular measurement, polygon measurement, curve measurement, ellipse measurement, arc measurement, point to point measurement, point to line measurement, parallel line measurement, angle measurement.

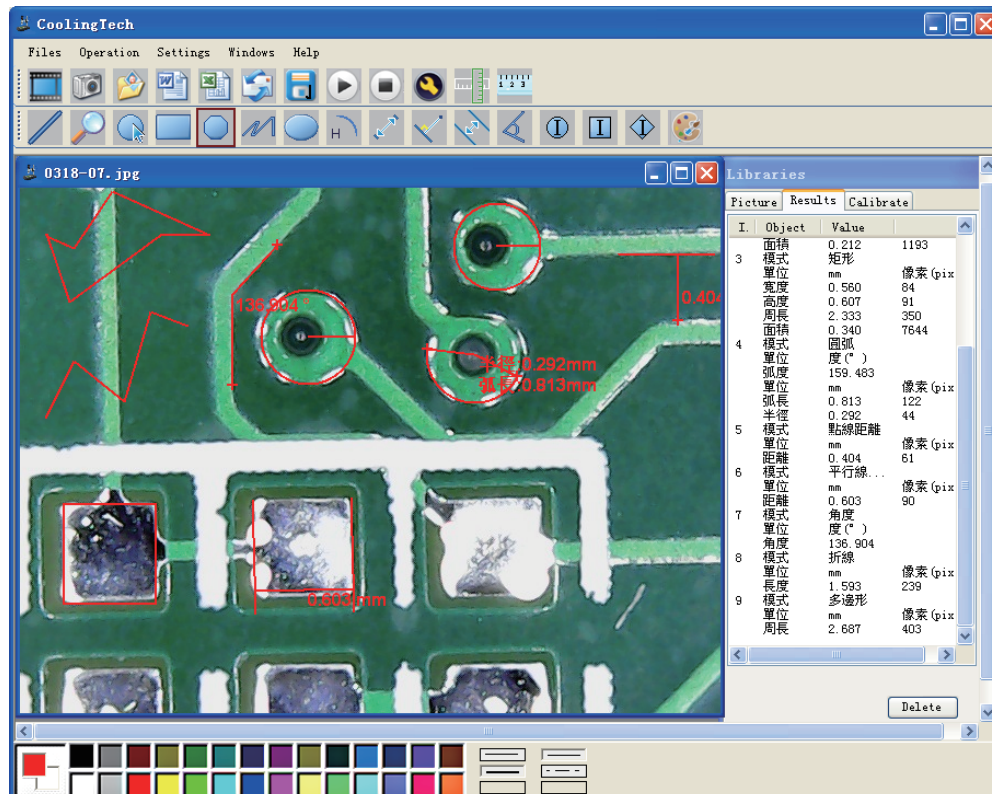

For example 1
Line measurement:

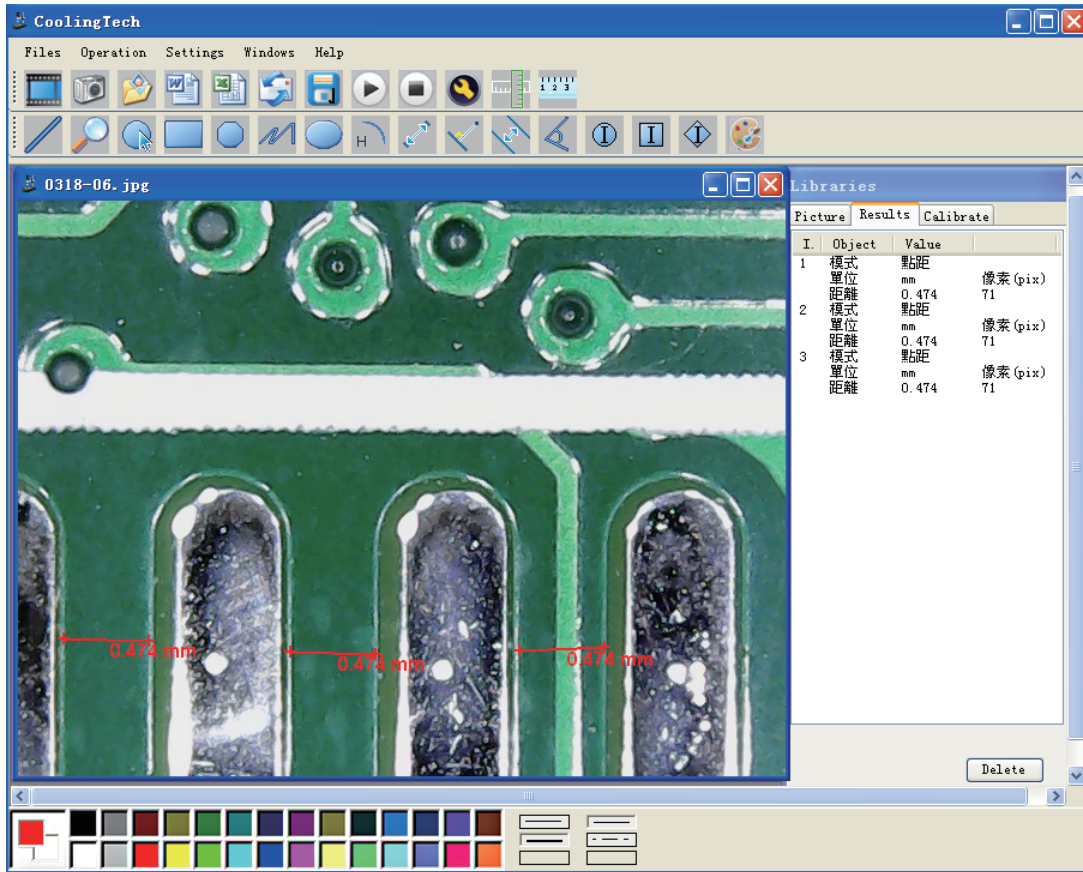

For example 2
Rectangular measurements:

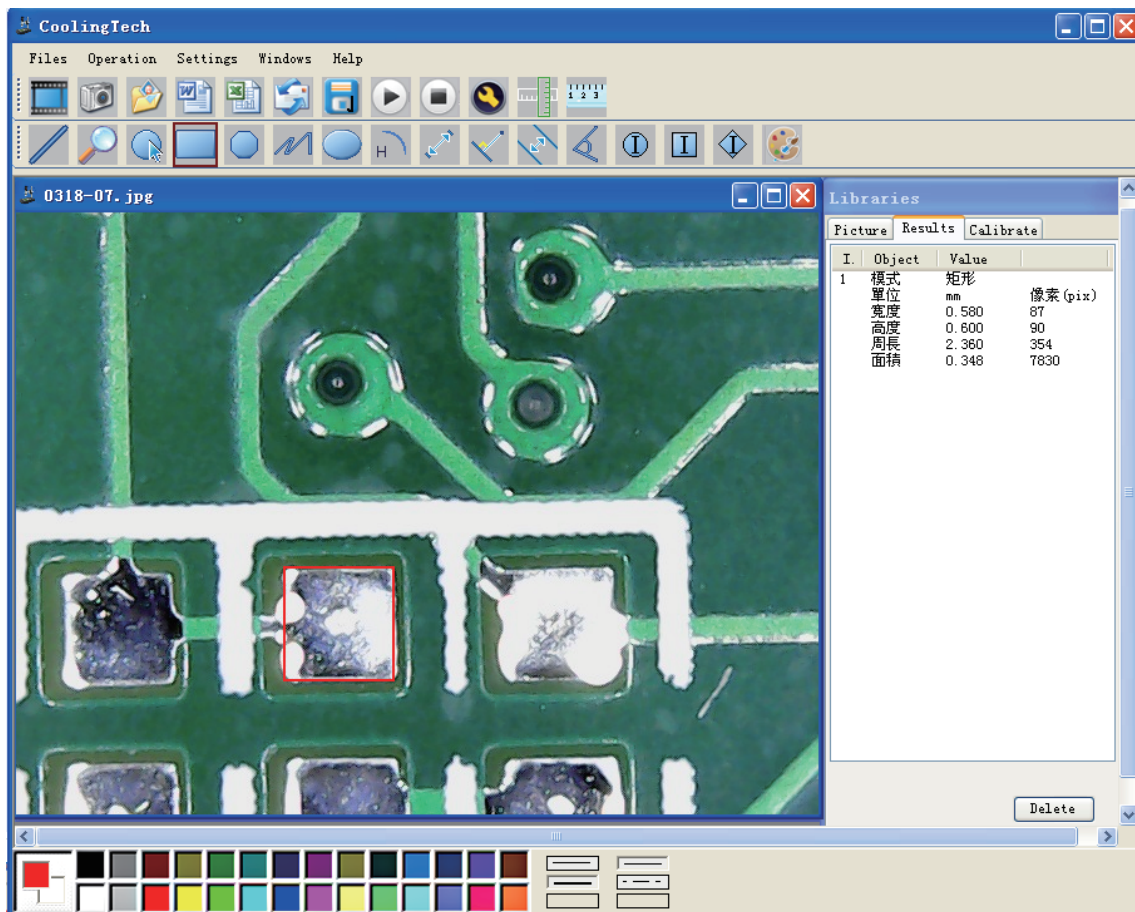

For example3
Ring measurements:

The screenshot shows the CoolingTech software interface. The main window displays an image of a green material with several circular features. One of these features is highlighted with a red circle, and a red line indicates a measurement across its diameter. The software's menu bar includes 'Files', 'Operation', 'Settings', 'Windows', and 'Help'. The toolbar contains various icons for file operations, navigation, and measurement tools. The 'Libraries' panel on the right shows a table of measurement results for two objects.

I.	Object	Value	
1	模式	圓	
	單位	mm	像素 (pix)
	半徑	0.667	100
	直徑	1.333	200
	周長	4.189	628
	面積	1.396	7853
2	模式	圓	
	單位	mm	像素 (pix)
	半徑	0.000	0
	直徑	0.000	0
	周長	0.000	0
	面積	0.000	0